




VEET 2.1 Quick Reference Guide

This guide is an overview. Please see the VEET Manual for detailed instructions. Software and documents referenced in this guide are available for download from www.projectveet.com

Last Updated: 07/23/2025


1

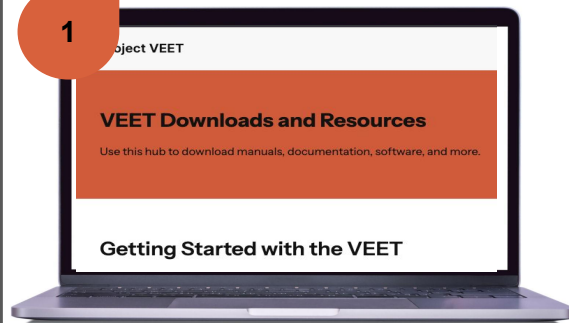

2



3

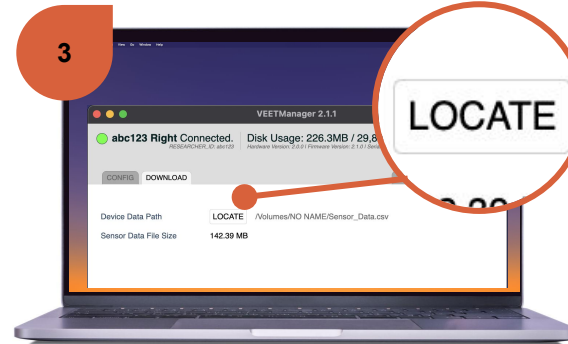

4



5



Setting up the VEET before handing the device to study participant

- Download and install the most recent version of the **VEETManager** from projectveet.com/downloads/
- Charge each VEET temple arm on wall power for at least 5 min.
- Connect one VEET temple arm to your computer. To avoid device conflicts, connect **only one** temple arm at a time.
- Open the **VEETManager** and upgrade to the latest firmware if prompted.
- Click the Download tab in **VEETManager** and Locate the VEET hard drive.
- Delete previous Sensor Data and Log Files.
- Adjust Participant / Researcher ID, Time Zone Offset, and sensor logging intervals on the **VEETManager** Configuration tab.
- Unplug and repeat the process with the opposite temple arm.

Fitting the VEET on study participant

- The VEET is compatible **only** with a few, specific glasses frame models.
- Please refer to the list of VEET-compatible glasses frames and order the most appropriate for your study participant.
- Replace the glasses frame temple arms with the VEET temple arms by removing and reusing the screws.
- Read the **Adjustable VEET 2.1 Optician's Guide** for instructions on how to assemble the Adjustable VEET, which comes with four pairs of temple tips and four pairs of hinges.


6




7




8

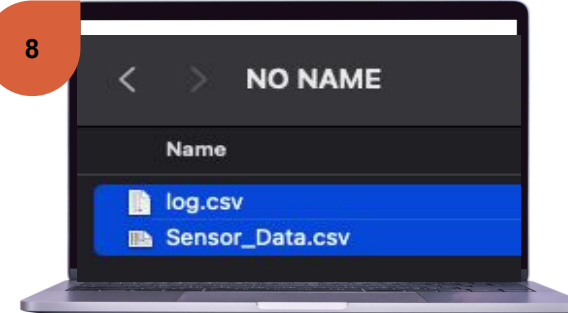

9

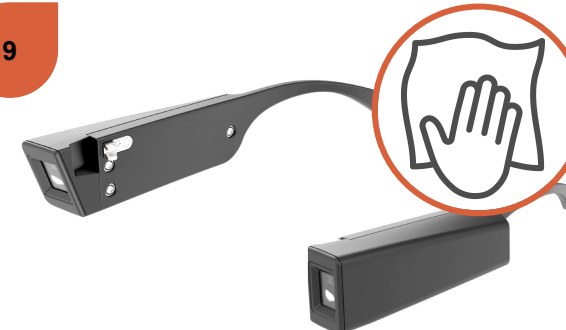

10

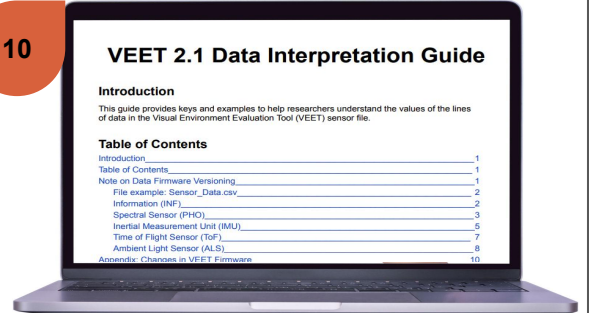

Unloading the data and preparing the VEET for reuse

- Before** handing VEET to study participants:
- Fully charge each temple arm for 3 hrs.
 - Connect the VEET to your computer and open **VEETManager** to ensure clock on the device is synchronized.
 - Instruct participants to charge VEET nightly on wall power. In default recording mode, VEET battery lasts 1.3 days.
 - Customize and provide **Research Participant Instructions (Template)**.
 - Take note of the date and time you provide VEET to study participant.
- After** retrieving VEET from study participants:
- Take note of the date and time you receive the VEET from the study participant.
 - Charge both VEET temple arms on wall power for at least 5 minutes.
 - Open the VEETManager and Locate the VEET hard drive.
 - Connect one VEET temple arm to your computer. To avoid device conflicts, connect **only one** temple arm at a time.
 - Copy the Sensor Data and Log Files to your computer.
 - Rename the files using helpful naming conventions (e.g. participant ID and "R" or "L" to indicate right / left temple arm).

- To prepare the device for reuse, delete any Sensor Data or Log files remaining on the VEET.
- Inspect the device for any physical damage.
- Wipe all surfaces with an isopropyl alcohol wipe.

Processing the VEET Data

- Validate your data by using the **VEET 2.1 Data Processing Guide**.
- Analyze your data by using the **VEET 2.1 Data Interpretation Guide**.
- Both guides are available for download from projectveet.com/downloads/